

VNMRJ Customization for APcheks

Place the file APchecks_good.tar in /home/username/vnmrsys directory and un-tar it. Files are copied into psglib, parlib, maclib, etc... All the file names in this pack start with AP so it is easy to list them for viewing purposes.

Issue the command *seqgen AP*.c* and this will compile the new pulse sequences.

cd to vnmrsys/templates/vnmrj/interface and issue *cp APMainMenuExpts.xml MainMenuExpts.xml*. This updates the pull-down Experiments menu in VNMRJ and shows the line AP checks of BP.

Currently Apchecks_good.tar is /home/vnmr2/vnmrsys and /home/vnmr3/vnmrsys directories in all the spectrometers.